

# Short Instructions *Seraline*<sup>®</sup>scan

Version: 5.0.1



1



prepare  
or edit  
Template

- Choose test
- ID: strip number
- Lab-ID: laboratory identification number
- Name of patient
- ↓
- Print template
- Lot.No: free selectable

2



- Fix developed strips on template

Strip No.	Controls	Serum Sample	Name	dsDNA	Nucleosomes	Sin	PO	Histones	U1 snRNP	RuSSA-A(500k)	RuSSA-A(250k)	LuSSB	Scl-70	CENP-B	Scl-1
1															
2															
3															
4															
5															
18															
19															
20															

3



new Scan

- Scan template
- Scanner glass should be tidy

4



Print Results

- Print results
- Table or single report